

Report

Capture Image

Export

Report


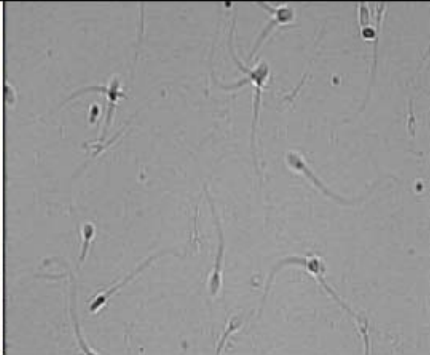
← BACK

After selecting either a series of records or one record, a report can be viewed, printed and/or exported to another location on the computer.

To run a **Stallion Semen Analysis Report**:

- Select the desired record by clicking in the record selection column.
- Click the REPORT button and the Stallion Report below will be displayed (the report will display the stallion picture and semen if they are attached to the record):

The Stallion name, picture and comments can be added to the report by going to: **Set-Up/Data Settings/Stallion Settings** (please refer to the Set-Up section of this User Guide)

STALLION SEMEN ANALYSIS REPORT			
Testing Facility			
Facility Name	Superior Stallions Stud		
Address	112454 Equine Drive	City	Cheyenne
State / Province	Wyoming	ZIP / Country Code	82001
Phone	+ 001 877-9090	Fax	+ 001 877-9091
E-Mail	mr_ed@superiorstud.com	Website	www.superiorstud.com
Stallion Information			
Name	Mr. Ed	ID/ Studbook #	2
Sample Processing			
Date	5/2/2008	Time	11:55
Sample #	21	Semen Volume [ml]	95.0
		Sample Type	Raw
TEST RESULTS			
Sperm Conc. [M/ml]	150.5	MSC [M/ml]	98.9
Motility [%]	65.5	PMSC [M/ml]	95.7
Prog. Motility [%]	63.6	Velocity [mic/sec]	89
Morph. [%]	69.4		
Totals per Semen Volume / Other Parameters			
# Sperm [Bil]	14.30	# Prog. Motile Sperm [Bil]	9.09
# Motile Sperm [Bil]	9.37	TNB	6499.21
Dosing Set-up			
Dosing Method	Prog. Motile Sperm	Maximum Dose Vol. [ml]	70.0
Minimum # Cells/Dose [M]	700		
Dosing Results			
Extender Volume [ml]	741	Dose Volume [ml]	69.6
# Doses	12	Total Volume [ml]	836
		# Sperm/Dose [M]	758
Comments			
Beautiful TV horse. Was on TV series: Mr. Ed - talking horse. Semen per straw \$675			
			

To run an SQA-Ve Test Report (with more than one test):

- Select the desired records by either clicking in the record selection column of the TABLE /setting a date range or selecting ALL.
- Click the REPORT button and the Report below will be displayed

SQA-Ve Test Report COOLED SEMEN

From: 4/27/2008 To: 4/28/2008

This sample SQA-Ve Test Report is for COOLED SEMEN by date range. Reports for all of the types of semen can be run based on the user selection.

SAMPLE DATA						TEST RESULTS							
Date	Time	Stallion		Sample #	Semen Volume [ml]	Sperm Conc. [M/ml]	Motility [%]	Prog. Motility [%]	MSC [M/ml]	PMSC [M/ml]	Velocity [mic/sec]	Totals per Semen Volume	
		ID/ Studbook #	Name									# Motile Sperm [Bil]	# Prog. Motile Sperm [Bil]
4/28/2008	03:34	8	Comanche	69	15.0	66.6	53.0	48.6	35.3	32.4	61	0.53	0.49
4/28/2008	03:29	25	Abaccus	68	17.0	51.5	84.1	32.7	43.3	16.8	86	0.74	0.29
4/28/2008	03:24	26	Bambi	67	20.0	83.4	76.8	67.4	64.1	56.2	89	1.28	1.12
4/28/2008	03:19	27	Lass	66	35.0	52.9	75.2	31.4	39.8	16.6	82	1.39	0.58
4/28/2008	03:14	28	Quark	65	30.0	49.9	70.6	29.7	35.2	14.8	69	1.06	0.44
4/28/2008	03:09	29	Magnum	64	20.0	54.6	72.7	31.0	39.7	16.9	83	0.79	0.34
4/28/2008	03:04	7	Trigger	54	16.0		<= 30					0.14	0.13
4/28/2008	02:59	46	Tex	53	20.0	49.9	42.9	31.6	21.4	15.8	55	0.43	0.32
4/28/2008	02:54	9	Jack	52	15.0	53.2	52.3	34.8	27.8	18.5	63	0.42	0.28
4/28/2008	02:49	48	Onix	51	20.0	73.8	73.3	61.0	54.1	45.0	79	1.08	0.90
4/28/2008	02:44	30	Mac	45	20.0		<= 30					0.06	0.05
4/28/2008	02:39	24	Bar Mount	44	22.0		<= 30					0.04	0.04
4/28/2008	02:34	42	Eagle Eye	43	15.0	55.9	40.9	23.5	22.9	13.1	53	0.34	0.20
4/28/2008	02:29	43	Oasis	28	25.0	50.9	36.7	24.1	18.7	12.3	50	0.47	0.31
4/28/2008	02:24	44	King	34	20.0	50.8	75.3	30.7	38.3	15.6	75	0.77	0.31
4/28/2008	02:19	45	Fabio	28	30.0	52.7	50.5	34.1	26.6	18.0	61	0.80	0.54
4/28/2008	02:09	39	Eddy	27	18.0	59.7	53.5	25.5	31.9	15.2	64	0.57	0.27
4/28/2008	02:04	40	Dancer	26	15.0	63.3	68.3	41.6	43.2	26.3	79	0.65	0.39
4/28/2008	01:59	41	Ted	22	17.0	55.1	50.3	25.8	27.7	14.2	61	0.47	0.24

This document was printed on 5/11/2009 17:08 from the E-Sperm system ver. 1.00 by A-Tech



- Click the Printer icon to print the report
- Click the Export button to save the data in Excel format (Excel software required)
- Use the ZOOM to minimize/maximize the view
- Use the page bar at the bottom of the report to move between pages
- Click the "X" in the upper right hand corner to exit